

Repair Enquiry Form for STable (©) Vibration Isolation Tables and Quiet Air Compressors

Date of request:

Contact person:

Mobile phone number of contact person:

Email address of contact person:

Location of the faulty equipment:

Type of repair action (indicate one option please):

Warranty repair

Extended Warranty repair

After warranty (invoiced) repair

What equipment went wrong? (indicate one option please):

STable (©) Vibration Isolation Table

Quiet Air Compressor

Name of Customer:

Date of shipment:

Serial number:

Table size:

Is the compressor working continuously? (you can hear it and there is a LED indicator on its front):

Which leg of the table does not keep its optimal working position? :

What is the position of the faulty working cylinder? (indicate one option please):

It is too high

It is too low

Can you see the mains switch of the Quiet Air Compressor lighting red, if the Compressor is switched on? :

What pressure can you see on the pressure meter of the Quiet Air Compressor, if the Compressor is switched on? :

Can you hear air leak noises? :

If you are able to locate the place of air leakage, where it is? :

Other remark or note:

If the compressor is warm, has a smell or generates a smoke, switch it off immediately!

If the compressor works continuously, please switch it off after the pressure check, not to override the 1000 hours of lifespan.